


SR8-Plus High Stand

Operation Manual

73-111-802

REV. A
DATE: 1998-12

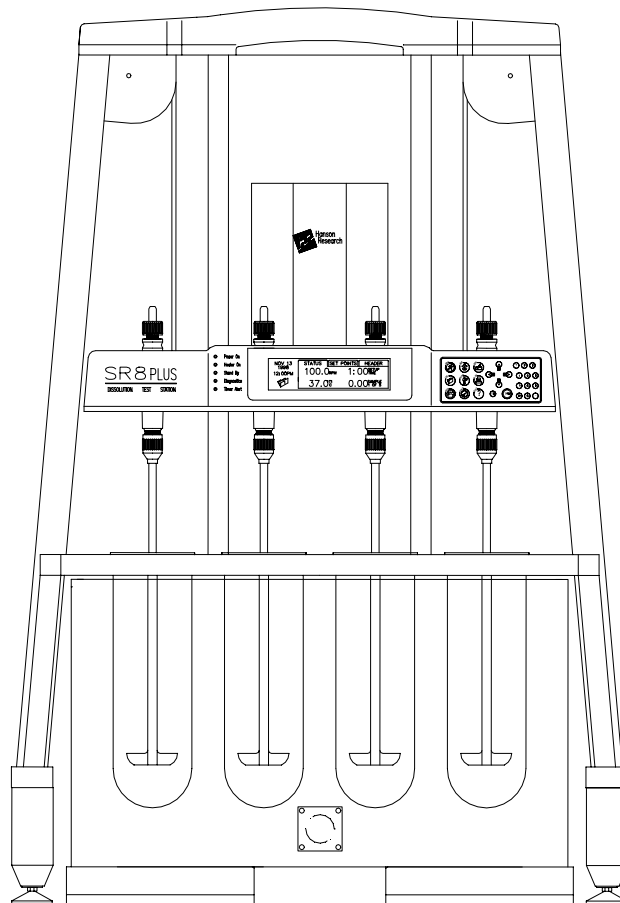

Hanson Research Corporation
9810 Variel Avenue • Chatsworth, CA 91311 USA
(800) 821-8165 • (818) 882-7266 • FAX (818) 882-9470
www.hansonresearch.com

©2004 Hanson Research Corporation.

Section One - Safety Considerations

General

1. Review all safety and environmental precautions pertaining to any chemicals, which are to be used in conjunction with this product.
2. Always wear eye protection.

Section Two - Introduction

Product Overview

The SR8-Plus High Stand is used for both 1 and 2 liter vessels. Legs are added to the unit to allow the taller bath to be placed under the base plate.

Section Three - Specifications

Follow the specifications listed in the main SR8-Plus operation manual with these exceptions.

Weight:

Dry:	68 kg (150 lbs.) – 115V
	70 kg (155 lbs.) – 230V
With liquid:	150 kg (283 lbs.) – 115V
	152 kg (288 lbs.) – 230V

Size:

Height: 112 cm (44 inches)

Bath Capacity:

With eight 1 liter flasks 61 liters (16 gallons)
With eight 2 liter flasks 53 liters (14 gallons)

Section Four - Installation

Legs:

Caution the unit is very heavy. Use at least two people to install legs.

1. Remove bath, bath support and control from back of unit.
2. The unit must be tipped to one side to install legs. Tip to one side and remove black feet.
3. Install these feet on the angled side of the new legs.
4. Install set screw into hole where feet were, turning screw in a few turns.
5. Screw leg onto setscrew so that the bottom of the stand comes into contact with the leg. Grab leg with two hands and tighten as much as you can, only tighten legs hand tight.
6. Repeat on other side.

Bath:

1. Follow instructions in SR8-Plus main manual. Slide bath all the way in, make sure to keep bath centered on the bath support.
2. Level unit per instructions in main manual.
3. Slide white plastic channels up to contact the bottom of the base, this will help reduce bath water evaporation.

Follow directions in the main SR8-Plus manual for all other installation and operation instructions.